Leica EG1120 – Instrucciones de uso

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Tecla RUN/STOP (CON/DES)

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RUN/STOP

MODE

Read the instructions for use carefully prior to operating the instrument, paying particular attention to the safety precautions!

2. **Connect** the power cable.
3. Turn on the mains switch (rocker switch with integrated green lamp). The **green POWER LED** on the control panel will illuminate. The POWER LED also indicates the standby condition.
4. Push the **RUN/STOP** button on the control panel in the standby to display the actual temperature of paraffin tank.
5. Turn the turning knob TANK to select the required set temperature (ambient to 70°C) of the paraffin tank.

There are two turning knobs on the control panel. Turn the **TANK** knob to select the set value of the paraffin tank (ambient to 70°C). Turn the **PLATE** knob to select the set value of the hot plate (ambient to 70°C).

The set values will be maintained, even if the instrument is switched off (RUN/STOP or mains switch), disconnected from mains or in the event of a power failure, provided the turning knob adjustment remains unchanged.

6. The display always reads the actual temperature of the paraffin tank.

Push the **MODE** button to change from actual paraffin tank temperature to actual hot plate temperature. The temperature of the hot plate will be indicated until the **MODE** button is released.

7. **To switch off** push the **RUN/STOP** button (standby). In daily routine it is not necessary to switch the instrument on and off with the mains switch.

8. Cleaning:

Owing to its smooth surfaces, the HistoTap is easy to clean. All surface materials are resistant to detergents commonly used in laboratories. The instrument must be switched off and cool for cleaning.

Organic solvents such as xylene must not be used for cleaning!

Safety precautions

Read the instructions for use carefully prior to operating the instrument and make yourself familiar with technical details of this instrument.

- Make sure that the power supply conditions in the laboratory correspond to the power rating indicated on the name plate of the instrument before connecting it to mains.
- The surface of the hot plate can be very hot during operation.

The instrument is designed for use with paraffin only!

Risk of burning!

Combustible or inflammable substances must not be stored near the instrument!

- The instrument may only be operated when filled. Minimum filling level: 1 cm
- For any repair and maintenance work, the instrument must be switched off and disconnected from mains.

Technical specifications subject to change.

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Leica Biosystems Nussloch GmbH