Leica EG1130 – Instrucciones de uso

Tecla RUN/STOP
(CON/DES)

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Tensión nominal:	100V, 60 Hz 120V, 60 Hz 230V, 50 Hz 240V, 50 Hz
Medidas:	460 x 235 x 625 mm
Arbeitsfläche:	290 x 460 mm
Gama de temperaturas de funcionamiento:	-5 bis -15 °C
Humedad del aire relativa:	10-80 %, sin condensación
Temperatureinstellbereich:	-5 °C bis -15 °C

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Leica EG1130 – Instructions for use

Technical Data

Nominal voltage:	100V, 60 Hz 120V, 60 Hz 230V, 50 Hz 240V, 50 Hz
Dimensions:	18,11 x 9,25 x 24,61 inches
work Dimensions:	11,42 x 18,11 inches
Dry weight:	77,09 lbs
Temperat. setting from:	-5 to -15 °C
Relative humidity:	10-80 %, non-condensing
Temperature control range:	room temperature up to 75 °C

Do not use sharp-edged tools to remove paraffin from the surface of the cold plate. Please use a soft plastic spatula only!

Safety precautions

The instructions for use must be read carefully prior to operating the instruments.

- It must be ensured that the power supply conditions in the laboratory correspond to the data indicated on the label of your Leica EG1130 prior to connecting it to mains. The instrument should be connected to shockproof sockets only.
- The instrument must be set up at a 10 cm distance on either side to ensure perfect operation of the cooling system.
- Prior to repair or maintenance, the instrument must be switched off and disconnected from mains.

1. **Connect** the power cable.
2. **Turn on** the main switch on the back of the instrument. The green **POWER** indication on the control panel at the front will illuminate.
3. Push the **RUN/STOP** button on the control panel (standby). The display reads the actual value of the cold plate.
4. The required **work temperature** is selected with the **potentiometer control knob**. The set value will be maintained when the instrument is turned off (RUN/STOP button or main switch) or even when mains power fails or if the instrument is disconnected from mains provided that the control knob position remains unchanged.
5. The Leica EG1130 is **switched off** using the **RUN/STOP** button. It is not necessary, however, to turn it off with the main switch in daily routine operation.
6. **Cleaning:**
The instrument has smooth surfaces. The control panel is protected with a foil. This ensures smooth, easy-to-clean surfaces. All surface materials are resistant to detergents commonly used in laboratories. The instrument must be switched off prior to cleaning.

Technical specifications subject to change.

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Leica Biosystems Nussloch GmbH