


Spectral Detector SP[®]

Leica TCS SP5 II: The Broadband Confocal
Leading in Multispectral Scanning

Living up to Life

Leica

MICROSYSTEMS

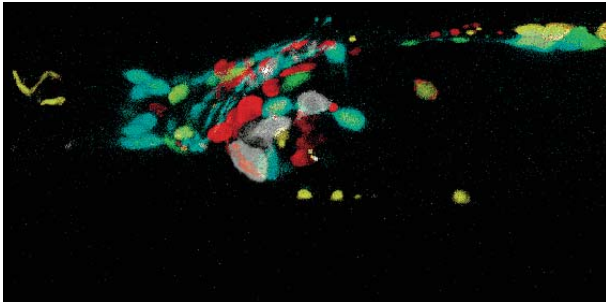

Caenorhabditis elegans, neurons.
5 channel image, overlay

- 1) Sensory and interneurons, CFP
- 2) Excitatory motoneurons, sensory neurons, YFP
- 3) GABAergic neurons, GFP
- 4) Glutamatergic interneurons, DsRed
- 5) Sensory neurons, DiD

Courtesy of Dr. Harald Hutter, Max Planck Institute
for Medical Research, Heidelberg, Germany

Multiparameter fluorescence





Fluorescence microscopy has become the major tool in modern biology. Nearly all disciplines in the biomedical field benefit from fluorescent staining. For structural analysis in morphology and anatomy, for kinetic measurements in physiology and biophysics, and for tracing gene-expression and dynamics in molecular biology – just to name the larger boundaries. Immunohistochemical fluorescence tags, fluorescence-in-situ hybridization, metabolite probes like Ca^{2+} or potential sensitive dyes and various specific stains for cell compartments and structural proteins fill the huge catalogues of fluorescence probes for current research tasks. And fashionable staining tools like fluorescent proteins and quantum dots flood the laboratories.

In addition to the great variety of dyes and probes there is another common trend in research: multiparameter fluorescence – application of multiple dyes to the same sample. This is an essential technique when interaction of proteins is the target. It is indispensable when communication between several individual cells is investigated and to find answers to questions in embryology – a classical example of identification of neuronal development by fluorescent proteins.

Colors Count

Past: cascade filtering – a nightmare for engineers and researchers



Of course, multi labeling is not a new invention. It was performed in the past by sequentially imaging the different channels. Or – if simultaneous recording was unavoidable – by cascading beam splitters and barrier filters. Whereas the first method is simply time consuming and prone to channel misalignment, the latter suffers from bulky filter magazines and still very limited flexibility. As a rule, the appropriate filter was simply not available – or the installation required a service call, too late for the already set up experiment and quite expensive, above all.

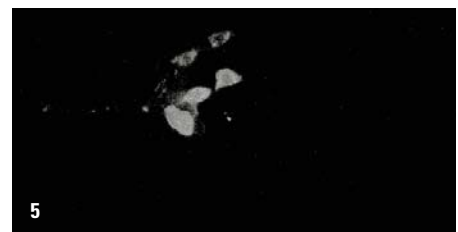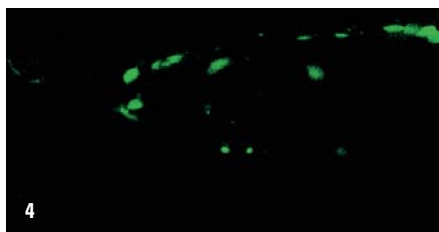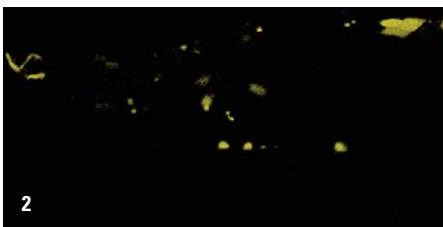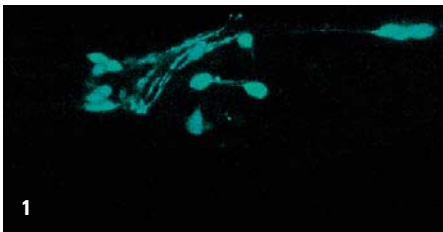

What is most unattractive with filter cascades is the low efficiency. Imagine constructing a cascade offering five channels. This system would at least have one channel where the light passes three different dichroic mirrors and a barrier filter. You can easily suspect that not too many photons would make their way through all these obstacles. As a result, the images are noisy – alternatively the sample bleaches, if you either average or increase the illumination intensity. Beside these efficiency problems, filters have a very strong limitation: they are not tunable. To adapt the system to various dye combinations, you are limited to the given characteristics of the filters and dichroic mirrors of your equipment. Although there might be combinations that “work”, they will seldom lead to optimal results due to loss in efficiency or unacceptable crosstalk – usually both.

Today: the Leica SP Detector

The Leica SP® detector solves all these problems. The light from the sample, which is a sum of multiple emission spectra, first passes through the pinhole. Here, the optical sectioning occurs. Then, by means of a prism, the colors are spread into a spectrum. On its way to the first detector, the light passes a slit photometer device consisting of two motorized barriers, which are controlled via software. The principle is simple but most effective: The device is steplessly adjustable: the barriers can be moved to any wavelength position. By means of this design, this single detector can make up any conventional filter specification in terms of center frequency and bandwidth.

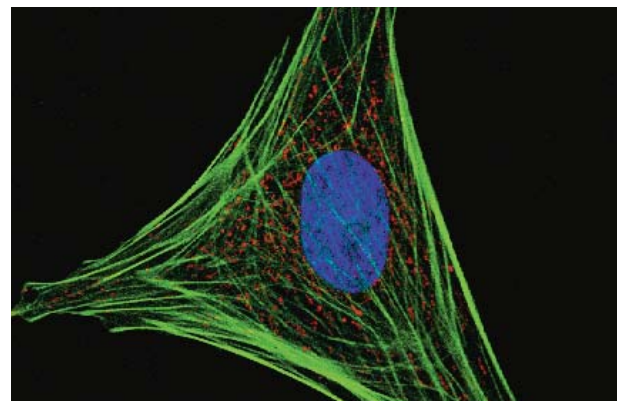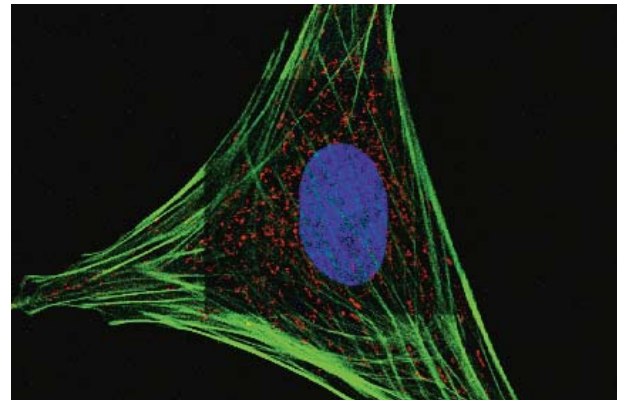

Sample protection by high sensitivity detection.


Top: high laser power and rigorous averaging is necessary in conventional detection systems in order to generate noise-free images.

Bottom: low illumination intensity and few averages will not damage this extremely bleach-sensitive sample when imaged with SP(R) detector.

An Ingenious Invention

However, this is not a multi-parameter fluorimeter. How can we simultaneously collect multiple segments from the spectrum? The clever bits are the barriers: they are not made of low-reflection material as it is the case for standard photometers. Here, the barriers are in fact mirrors. So the residing parts of the spectrum on both sides of the band collected at sensor 1 are directed to sensors 2 and 3 with again motorized, reflective barriers – ready for the next sequence.

This is how the SP5 II detector works. Five channels simultaneously. Full efficiency. No filters and dichroics. Just some high-efficiency coated mirrors. And all bands unrestrictedly tunable. Whether you need to cover the full spectrum for full efficiency or just small specific bands to the detector for unsurpassed crosstalk reduction, the decision is yours! (Software for unmixing is, of course, available anyway.)

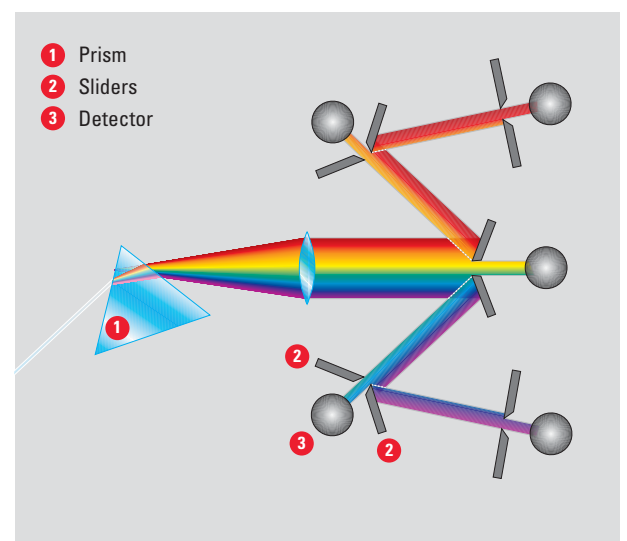

Better images in less time

Your benefit from Leica's SP® detector concept: maximum efficiency and full tunability which guarantees you best imaging quality and least crosstalk in multiparameter fluorescence. And as a side-effect: the SP® detector allows recording of lambda scans – spectral image stacks to analyze the in-situ emission characteristics of the dyes employed in your sample. There have been several attempts to top the versatility of this concept, but mainly targeting the issue of spectral imaging. Some approaches work with optical grids instead of prisms – by far too inefficient for imaging. There are attempts to overcome the grid problem, only by introducing even more optical elements, consequently eating photons. And cumbersome serial-filter systems can do basic spectral scanning, but at very limited accuracy.

FRET

After all, the real benefit of the SP® detector is not lambda scanning, but efficiently collecting specific emissions from different dyes in the sample. Most applications would not even need lambda scanning. For example, if you want to monitor FRET it is sufficient to monitor just two channels, since the changes occur spectrally insensitively and only one species increases, whereas the other decreases. All you need is two channels – however, with the correct band settings.

All Those Benefits

How the SP detector fits the future

Confocal microscopes – even when equipped with an SP® detector – still suffer from two main obstacles. Firstly, there is still the primary beam splitter: the element which splits the excitation for illumination and the emission for detection. Classically, this is again a device which is not tunable and not very efficient, either. The Leica AOBS® overcomes these restrictions. The AOBS®, an acousto optical beam splitter, is a tunable element at very high transmission, which switches within a matter of microseconds to allow you to appropriately channel the illumination and emission bands.



- **Spectral Detector, five confocal channels**
- True confocal point-scanning – real optical sectioning
- High dynamic sensors
- Low dark current PMTs
- High efficiency prism spectrometer
- Loss-free band selection
- Free tunable emission bands
- Any dye adaptable
- Five bands simultaneously
- Dye separation: sufficient for five channels simultaneously

Optionally, Leica Microsystems offers two fast sensors connectable to detection equipment for fluorescence lifetime imaging within the SP® detector. This combination allows FLIM measurements to be performed in any emission segment of the spectrum. A sequence of FLIM measurements at incremented wavelengths provides a new dimension in microscopy: fluorescence lifetime as a function of emission color “SP FLIM”. A new area, which is still waiting for exploration. Completely new applications will be possible. More differentiated insights in FRET processes by simultaneous detection of acceptor and donor, and other quenching effects are now available for research. In addition, SP-FLIM has better photon counting efficiency than filter-based concepts.

Spectral FLIM – a New Dimension

Fluorescence Lifetime Imaging

To further increase the range of applications, it is possible to install two different sources for FLIM excitation in the same system: either a pulsed Ti:sapphire laser, which is the standard source for multiphoton excitation, or a pulsed 405 nm laser diode.

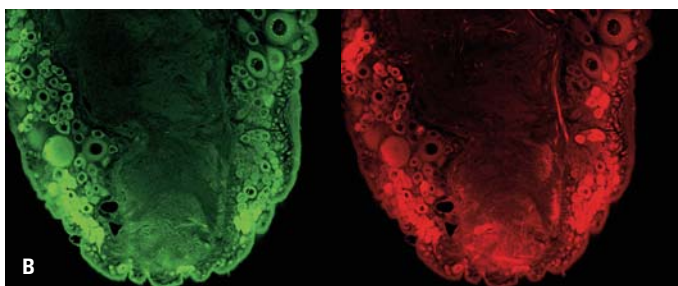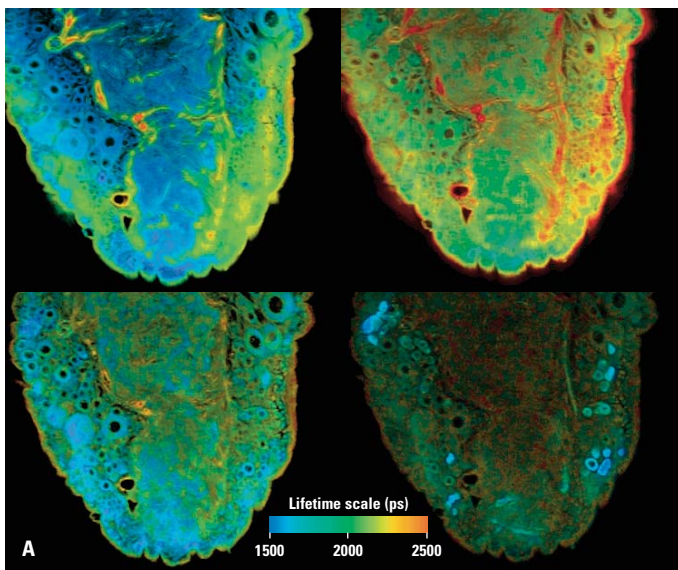

Schistocerca gregaria, nervous system, double staining Cy2, Cy3.

A) SP-FLIM images

left: emission 500-550 nm, right: emission 575-647 nm

top: D-FLIM, excitation 405 nm, bottom: MP-FLIM, excitation 780 nm

B) Intensity images for comparison

left: excitation 488 nm, emission 500-550 nm

right: excitation 561 nm, emission 575-647 nm

“With the user, for the user”

Leica Microsystems

Leica Microsystems operates globally in four divisions, where we rank with the market leaders.

● Life Science Division

The Leica Microsystems Life Science Division supports the imaging needs of the scientific community with advanced innovation and technical expertise for the visualization, measurement, and analysis of microstructures. Our strong focus on understanding scientific applications puts Leica Microsystems' customers at the leading edge of science.

● Industry Division

The Leica Microsystems Industry Division's focus is to support customers' pursuit of the highest quality end result. Leica Microsystems provide the best and most innovative imaging systems to see, measure, and analyze the microstructures in routine and research industrial applications, materials science, quality control, forensic science investigation, and educational applications.

● Biosystems Division

The Leica Microsystems Biosystems Division brings histopathology labs and researchers the highest-quality, most comprehensive product range. From patient to pathologist, the range includes the ideal product for each histology step and high-productivity workflow solutions for the entire lab. With complete histology systems featuring innovative automation and Novocastra™ reagents, Leica Microsystems creates better patient care through rapid turnaround, diagnostic confidence, and close customer collaboration.

● Surgical Division

The Leica Microsystems Surgical Division's focus is to partner with and support surgeons and their care of patients with the highest-quality, most innovative surgical microscope technology today and into the future.

The statement by Ernst Leitz in 1907, “with the user, for the user,” describes the fruitful collaboration with end users and driving force of innovation at Leica Microsystems. We have developed five brand values to live up to this tradition: Pioneering, High-end Quality, Team Spirit, Dedication to Science, and Continuous Improvement. For us, living up to these values means: **Living up to Life.**

Active worldwide

Australia:	North Ryde	Tel. +61 2 8870 3500	Fax +61 2 9878 1055
Austria:	Vienna	Tel. +43 1 486 80 50 0	Fax +43 1 486 80 50 30
Belgium:	Groot Bijgaarden	Tel. +32 2 790 98 50	Fax +32 2 790 98 68
Canada:	Richmond Hill/Ontario	Tel. +1 905 762 2000	Fax +1 905 762 8937
Denmark:	Herlev	Tel. +45 4454 0101	Fax +45 4454 0111
France:	Nanterre Cedex	Tel. +33 811 000 664	Fax +33 1 56 05 23 23
Germany:	Wetzlar	Tel. +49 64 41 29 40 00	Fax +49 64 41 29 41 55
Italy:	Milan	Tel. +39 02 574 861	Fax +39 02 574 03392
Japan:	Tokyo	Tel. +81 3 5421 2800	Fax +81 3 5421 2896
Korea:	Seoul	Tel. +82 2 514 65 43	Fax +82 2 514 65 48
Netherlands:	Rijswijk	Tel. +31 70 4132 100	Fax +31 70 4132 109
People's Rep. of China:	Hong Kong	Tel. +852 2564 6699	Fax +852 2564 4163
Portugal:	Lisbon	Tel. +351 21 388 9112	Fax +351 21 385 4668
Singapore		Tel. +65 6779 7823	Fax +65 6773 0628
Spain:	Barcelona	Tel. +34 93 494 95 30	Fax +34 93 494 95 32
Sweden:	Kista	Tel. +46 8 625 45 45	Fax +46 8 625 45 10
Switzerland:	Heerbrugg	Tel. +41 71 726 34 34	Fax +41 71 726 34 44
United Kingdom:	Milton Keynes	Tel. +44 1908 246 246	Fax +44 1908 609 992
USA:	Bannockburn/Illinois	Tel. +1 847 405 0123	Fax +1 847 405 0164

and representatives in more than 100 countries