# Leica HI 1210 – Instrucciones de uso

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RUN/STOP  
(CON/DES)

SET

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## Tastensymbole

RUN/STOP

Arrow buttons for temperature selection

SET

1. **Connect** the power cable.
2. **Turn on** the mains switch on the back of the instrument. The green **POWER** lamp on the control panel will illuminate.
3. Push the **RUN/STOP** button on the control panel to heat the Leica HI1210 up to the previously entered set temperature. The display reads the actual value. On reaching a temperature of approx. 44 °C, the yellow **DANGER/HOT** lamp illuminates. For reasons of safety, the yellow lamp illuminates continuously when a temperature of 44 °C is reached, not only during the heating periods.
4. The required temperature **is selected** using the **ARROW** buttons. At this stage, the indicated value is the set temperature. Once the appropriate temperature is displayed, the **ARROW** button should be released. The indicated set value will be memorised automatically. After about 2 seconds the display returns to actual temperature. The required temperature can be selected either step by step by repeatedly pushing the appropriate **ARROW** button, or continuously by holding the button down.
5. The set temperature can be verified by pushing the **SET** button. The set temperature is **maintained** even if the instrument is switched off (with the RUN/STOP button or mains switch), disconnected from mains, or in the event of a power failure.
6. **To switch off** push the **RUN/STOP** button. In daily routine operation, it is not necessary, however, to switch the instrument off with the mains switch.

## 7. Cleaning:

Lateral ventilation slits are avoided on the Leica HI 1210. The control panel is protected with a foil. This ensures smooth, easy-to-clean surfaces. All surface materials are resistant to detergents commonly used in laboratories. The instrument must be switched off and cool for cleaning.

**Organic solvents such as xylene must not be used for cleaning!**

## Safety precautions

The instructions for use must be read carefully prior to operating the instrument.

- It must be ensured that the power supply conditions in the laboratory correspond to the power rating indicated on the name plate of the instrument before connecting it to mains.
- The surface can be very hot during operation.

**Flammable liquids must not be used!**

**Risk of burning!**

**Combustible or flammable substances must not be stored near the instrument!**

- The instrument must be filled when operated. Minimum filling level: 1 cm
- Prior to repair or maintenance, the instrument must be switched off and disconnected from mains.

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Leica Biosystems Nussloch GmbH